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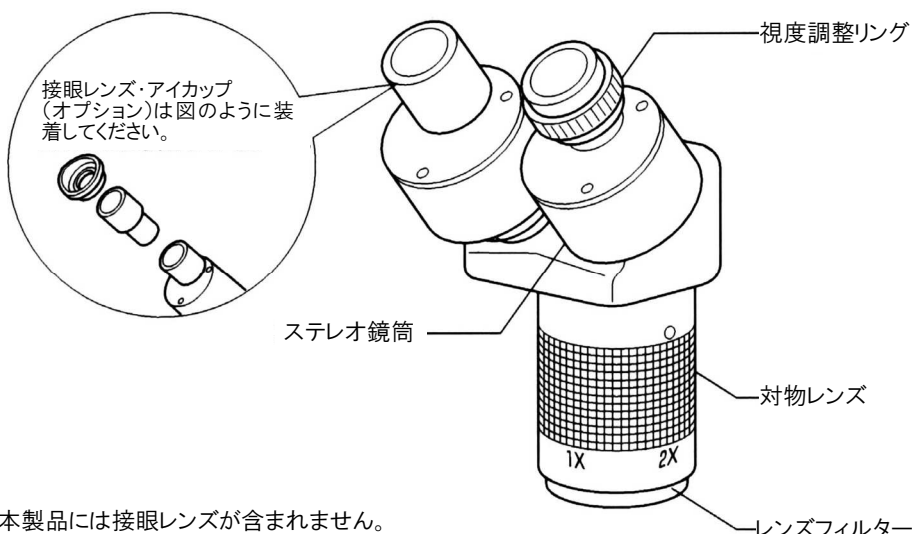
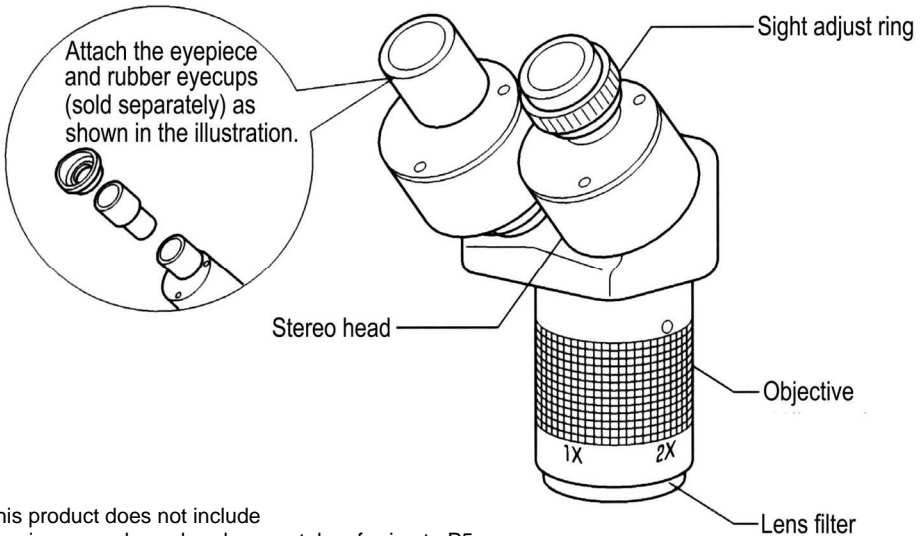
TEL(06)6567-3111 FAX(06)6562-0024


# L-501 STEREO HEAD SCOPE

Thank you for purchasing the HOZAN L-501 STEREO HEAD SCOPE. With proper care and handling, this fine instrument will provide years of trouble-free operation. Please read this entire instruction manual carefully before attempting to place this instrument in service. Please keep this instruction manual available for reference.

## Identification of parts

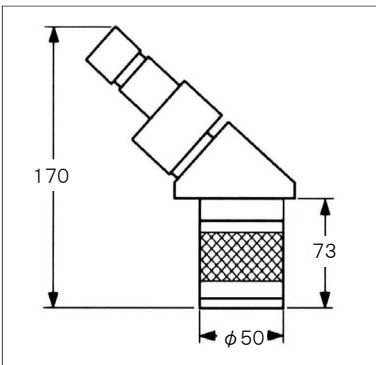\* This product does not include Eyepieces can be ordered separately referring to P5.



## Specifications

Objectives	1X、2X、 selection format
Eyesight adjustment	Monocular adjustment type, ±5Dptr
Working Distance	80mm
Focus adjustment	Head scope up/down type
Stereo head spec.	Binocular 45°tilted head scope
Pupil distance range	55-74mm · Independent left/right type

## Warning and caution symbols

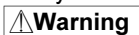These symbols are used throughout the instruction manual to alert the user to potential safety hazards as follows :

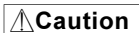 **Warning** ... Notice when incorrect handling could cause the user's death or serious injury.

 **Caution** ... Notice when incorrect handling could cause injury to the user or material damage.

Even if the instructions do not have  **Caution** mark, there are some possibilities for a serious situation. Follow the instructions.

## Precautions

The L-501 STEREO HEAD SCOPE is designed for precision inspection of PCB.  
Do not use for any other purpose.

### **Caution**

1. Use only genuine HOZAN component as the holder (e.g., L-509, L-522).
2. Do not unnecessarily remove the stereo head from the post. If it becomes necessary to remove the stereo head, hold it with both hands to prevent dropping the head. This could damage the head itself, base, or specimen.
3. Do not disassemble or modify the L-501 yourself, otherwise this could cause a malfunction.
4. Protect the unit against physical shock or vibration during transport. Otherwise, this could cause a malfunction.
5. Do not expose the unit to prolonged direct sunlight, high temperature, high humidity, or dust. Such exposure could cause a malfunction.

## Note

- Keep the lenses and filter clean from dust and fingerprint. Otherwise, this could cause an obscure image. (Ref. daily care on page 4.)

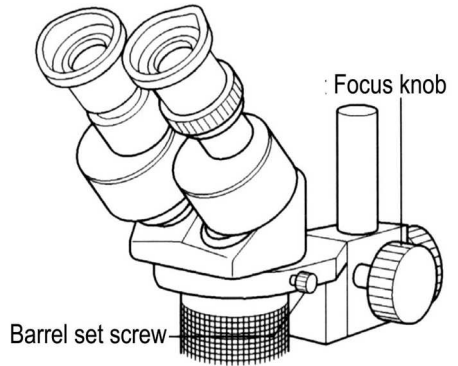
## Operation

- 1 Set up the base with an L-509 attached or an L-522 on a level and stable surface.
- 2 Insert the L-501 into the holder.


## Operation

- 3** Tighten the barrel set screw to secure the barrel to the holder.  
Attach the eyepieces (e.g.,L-511) and conversion lens.


- ※ The distance from the stage to the end of the objective is 80 mm. For the L-521, align the top edge of the L-509 holder with the matching mark on the post.

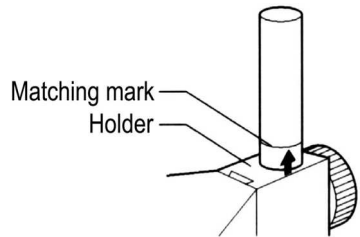- 4** Place the specimen on the stage.

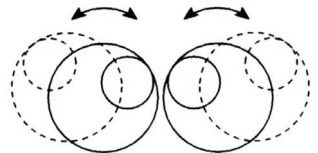
- 5** While looking through both eyepieces, turn the stereo barrels to view a single circle.

- 6** While looking only through the left eyepiece with your left eye, turn the focus knob to adjust the focus.

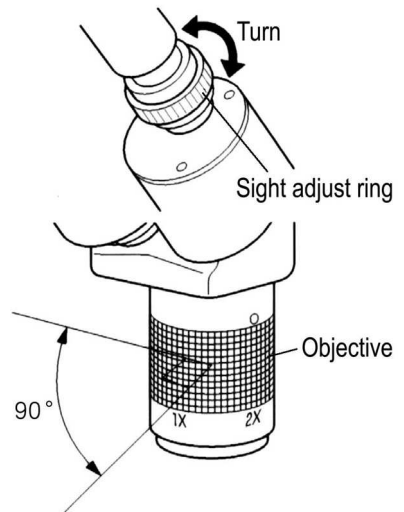- 7** Then, while looking only through the right eyepiece with your right eye, turn the sight adjust ring to adjust the sight.


- ※ The magnification can be switched between 1× and 2× by turning the objective. Since turnable angle of objective is 90°, if it cannot turn to one direction, turn it in the opposite.

## Operation
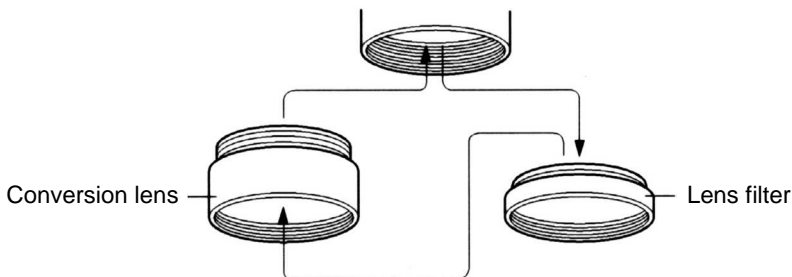
### ● Using the conversion lens

Rotate and remove the lens filter at the bottom and install the conversion lens onto the objective. Attach the lens filter at the bottom again.

※ Refer to the page 5 for details on the magnification, field of view and working distance when using the conversion lens.

## Daily care

### Lens

- Wipe away any dust gently with a soft cloth or a piece of gauze.
- Clean away fingerprints or oil stains with a cloth which is lightly soaked in denatured alcohol (such as ethyl alcohol or methyl alcohol).

※ Alcohol is extremely flammable, so handle it with care.

### Stereo head and plastic parts

- Wipe with a soft, dry cloth.  
Do not use any organic solvents such as thinner, alcohol or ether.
- When out of use for an extended period, store this unit in a dry dust-free place.

## Optional accessories

Use only the following optional accessories manufactured by HOZAN.

### Lighting systems

#### L-711 LED light

The specimen scarcely receives any temperature influence from this LED light, since the generation of heat is extremely insignificant by comparison with other illuminants. The light has few fluctuations of brightness and no flicker at all.

Various lighting systems can be used. Please refer to our catalog or web catalog.

# Optional accessories

## Photographic equipment

### L-835 USB camera

Specimens under a microscope are photographed (still and video) mounting on it through the L-845 Adapter kit. Still images can be edited using the provided software. Without lenses.

L-835

### L-845 Adapter kit

Use for connecting a camera which has C mount to a microscope.

L-845

## Eyepiece

Use either the L-511(10×), L-512(20×), L-513(15×).

These lenses can easily be removed just by pulling them out.

- Two per pack
- With eyecups

## Eyepiece with scale

- L-69(10×)

Use as replacement for monocular lenses only.

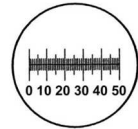
■ Magnification: 10x ■ Quantity: 1

\* Not usable where the total magnification is other than 10x.

\* An eye cup accessory is not provided.

L-69

Micrometer  
(5 mm by  
1/10 mm step)

## Conversion lens

Use either the L-57-0.5(0.5×), L-57-1.5(1.5×) or L-57-2.0(2.0×).

When installing, remove the accessory lens filter and screw on the conversion lens. Then re-attach the lens filter.

The magnification, field of view and working distance gotten when combining eyepieces, objectives and conversion lenses are follows.

### Magnification table

Standard item

Objective	1×			2×			1×			2×			1×			2×								
Conversion lens							L-57-0.5						L-57-1.5						L-57-2.0					
Eyepiece	Magnification ×	Field of view mmφ	Working distance mm	Magnification ×	Field of view mmφ	Working distance mm	Magnification ×	Field of view mmφ	Working distance mm	Magnification ×	Field of view mmφ	Working distance mm	Magnification ×	Field of view mmφ	Working distance mm	Magnification ×	Field of view mmφ	Working distance mm	Magnification ×	Field of view mmφ	Working distance mm	Magnification ×	Field of view mmφ	Working distance mm
L-511 (WF10×)	10	18		20	9		5	36		10	18		15	12		30	6		20	9		40	4.5	
L-513 (WF15×)	15	12	80	30	6	80	7.5	24	150	15	12	150	22.5	8	40	45	4	40	30	6	27	60	3	27
L-512 (WF20×)	20	9		40	4.5		10	18		20	9		30	6		60	3.6		40	4.5		80	2.2	

\*The "working distance" is the distance from the end of the filter to the subject.

**HOZAN TOOL INDUSTRIAL CO.,LTD.**  
1-2-12 Saiwaicho, Naniwa-ku, Osaka 556-0021, Japan  
Tel : 81-6-6567-3111 Fax : 81-6-6562-0024