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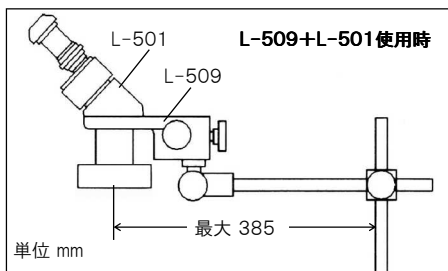
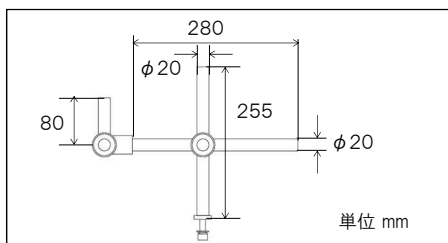
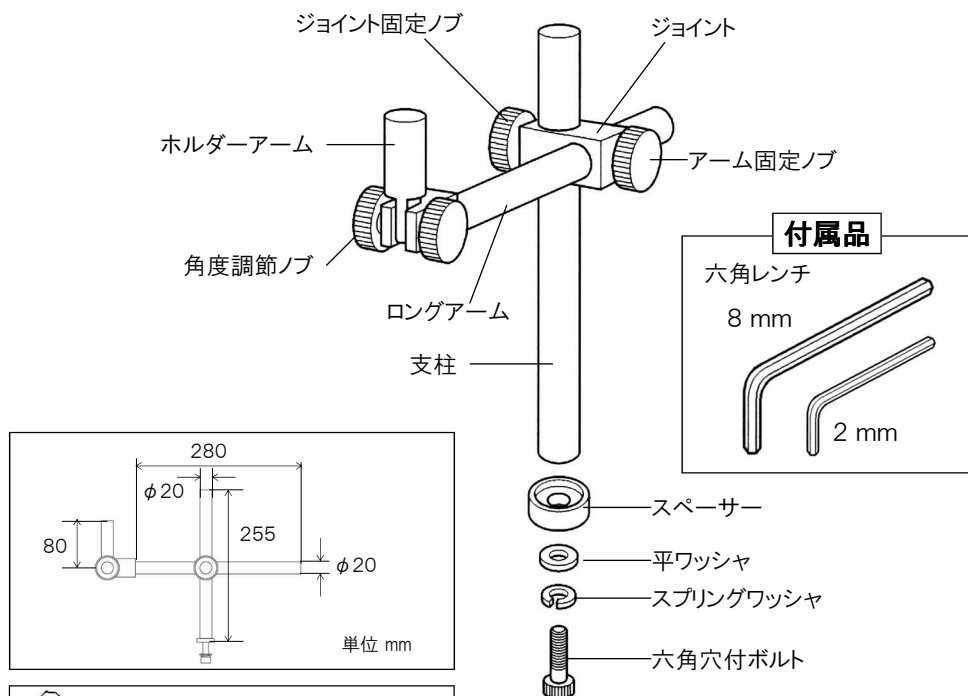
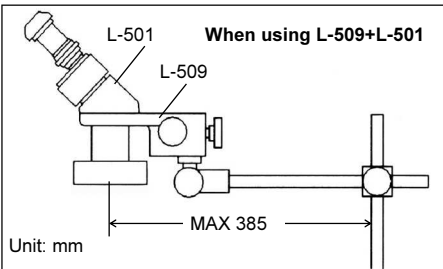
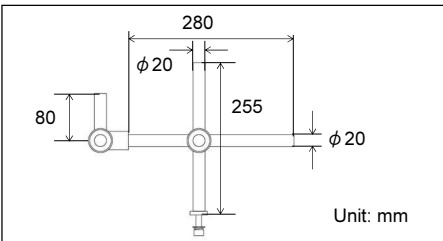
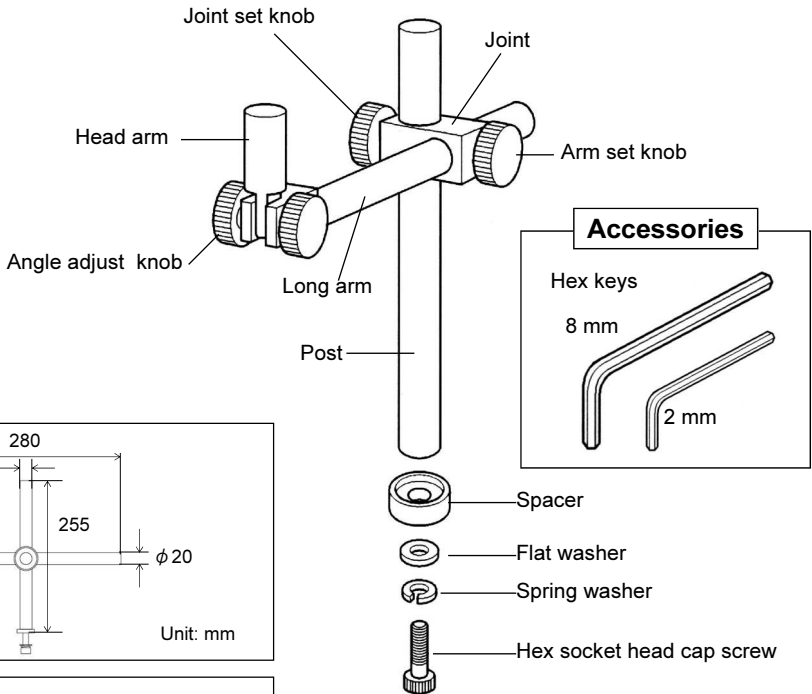
# L-523 LONG TYPE ARM

Thank you for purchasing the HOZAN L-523 LONG TYPE ARM. With proper care and handling, this instrument will provide years of trouble-free operation. Please read this entire instruction manual carefully before attempting to place this instrument in service. Please keep this instruction manual available for reference.

## Identification of parts and contents

## Specifications

Post diameter	20 mm φ
Post height	255 mm
Arm length	280 mm

## Warning and caution symbols

These symbols are used throughout the instruction manual to alert the user to potential safety hazards as follows :

**⚠ Warning** ... Notice when incorrect handling could cause the user's death or serious injury.

**⚠ Caution** ... Notice when incorrect handling could cause injury to the user or material damage.

Even if the instructions do not have **⚠ Caution** mark, there are some possibilities for a serious situation. Follow the instructions.

## Precautions

The L-523 is an arm for use only in installing the L-501 Stereo Head and such to the L-525 Large Table. Do not use for any other purpose.

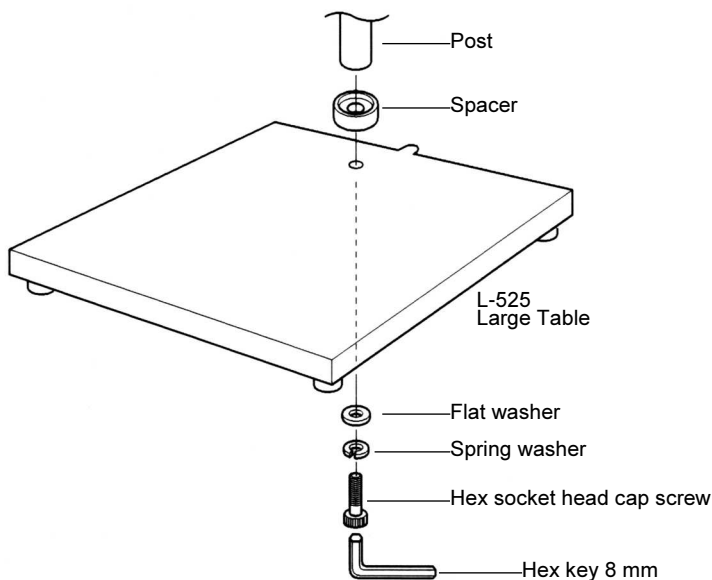
### ⚠ Caution

Be sure all parts are tightened securely. Otherwise, the microscope or the CCD camera could fall causing damage for them and the specimen.

## Operation

### Prepare the L-525 Large Table (sold separately) for installing the L-523.

- 1 Assemble the post, spacer, flat washer, spring washer, and hex socket head cap screw with the L-525 Large Table as shown in the illustration. Tighten the screw securely with the hex key 8 mm provided.

## Operation

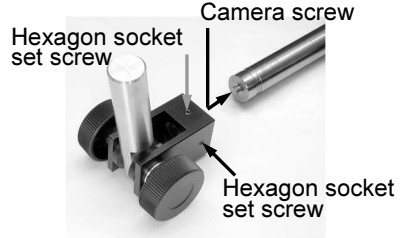
- 2 When use the ESD rubber mat, insert it into the post and place on the table, surely grounding. Refer to the instruction manual of the L-525 for detailed explanation.
  
- 3 Place the joint on the post, and then secure the joint set knob.
  
- 4 Insert the long arm into the joint. Then, secure the arm set knob after deciding the position back and forth.
  
- 5 Secure the angle adjust knob after deciding the angle of the holder arm as required.
  
- 6 Insert the optional head holder into the head arm, and secure the knob of the head holder.  
Select among L-509 Stereo Head Holder, L-809 Camera Platform with Slider, L-804 Micro-Drive Head Holder and such as a holder.  
Refer to HOZAN catalogue and our website for details.

# Operation


## Removing the holder arm section

A camera screw will appear when removing the holder arm section. A pan head or a camera in the market can be connected on this.


The holder arm section has been secured on the long arm with two hexagon socket set screws. To remove the holder arm section from the long arm, unscrew the hexagon socket set screws using the hex key 2 mm provided. Do not remove the screws; screw in them avoiding to lose them after removing the holder arm section.

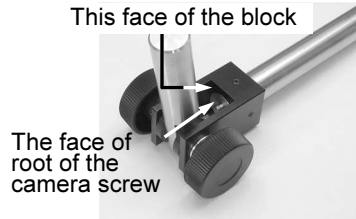To install the holder arm section again, secure with the hexagon socket set screws aligning the two faces explained in the illustration.

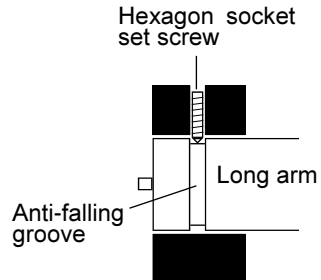Make sure that hexagon socket set screws are catching the anti-falling groove.

**HOZAN TOOL INDUSTRIAL CO.,LTD.**

1-2-12 Saiwaicho, Naniwa-ku, Osaka 556-0021, Japan  
Tel : +81-6-6567-3111 Fax : +81-6-6562-0024