L-850-2

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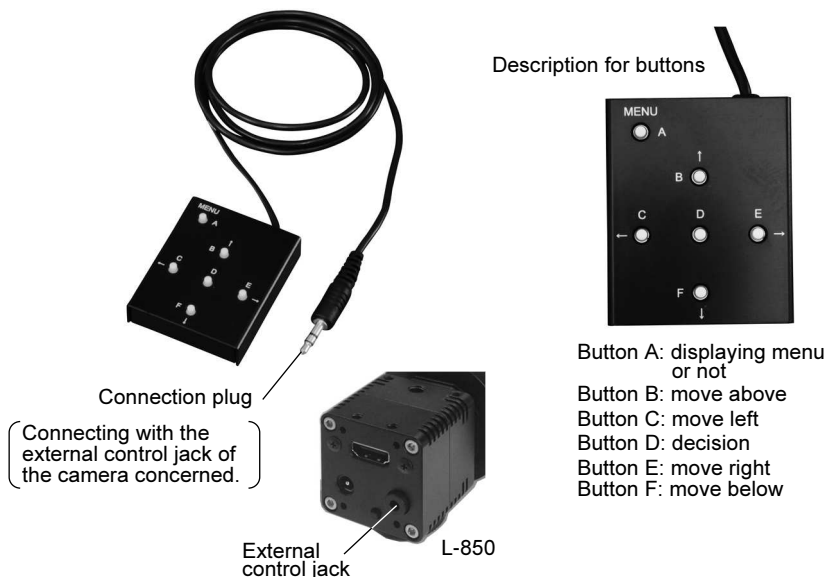
L-850-2 OSD CONTROLLER

Thank you for purchasing the HOZAN L-850-2 OSD CONTROLLER. With proper care and handling, this fine instrument will provide years of trouble-free operation. Please read this entire instruction manual carefully before attempting to place this instrument in service. Please keep this instruction manual available for reference.

Description for buttons

- Button A: displaying menu or not
- Button B: move above
- Button C: move left
- Button D: decision
- Button E: move right
- Button F: move below

Warning and caution symbols

These symbols are used throughout the instruction manual to alert the user to potential safety hazards as follows :

Warning ... Notice when incorrect handling could cause the user's death or serious injury.

Caution ... Notice when incorrect handling could cause injury to the user or material damage.

Even if the instructions do not have **Caution** mark, there are some possibilities for a serious situation. Follow the instructions.

Precautions

This device is for Hozan's L-850 FULL HD CAMERA. Do not use for any other purpose.

Caution

1. Do not disassemble or modify the device. This could cause malfunction.
2. Do not soak into water. Do not splash water.
3. Do not give a strong impact. This could damage the inner mechanisms.
4. Do not leave under the scorching sun or in an environment of high temperature.

Operation

Connect the connection plug to the external control jack of the target camera.

Setting window

Press the A button to display the setting menu on the computer screen.

PAGE 1 2 3 4 5 6

▶ AE MODE	NORMAL
LUMINANCE	00000
PRI GAIN	27.09dB
PRI SHUTTER	0000 1/1000s
USER GAIN	000 46.2dB
USER SHUTTER	000 1/1000s
GAMMA	0.45

Using the B/F button, move the cursor to each setting item. Then depress the E button. The letters of the item reverse negatively. Alter them using the B/F button. Return by the C button. The menu has six pages. Simply pass one after another by the B/F button.

Revised settings using this device can be saved in the camera. When saving new settings, former settings are not held. Please operate prudently not to delete the initial settings or the formerly saved settings referring "How to save" on P4.

Description for various settings

In [] are the initial settings.

PAGE 1 Setting brightness of the image

AE MODE Setting auto exposure control. **[NORMAL]**

NORMAL: Controlling exposure automatically.

LONG: Controlling exposure time over the frame rate. Enough brightness can be gotten even under an insufficient illumination.

USER MODE: Setting exposure manually. Adjust the values of USER GAIN and USER SHUTTER.

LUMINANCE: Setting luminance. **[4608]**

PRI GAIN: Setting gain. **[OFF]**

PRI SHUTTER: Setting exposing time. **[0000]**

USER GAIN: Setting gain when selecting USER MODE. **[000]**

USER SHUTTER: Setting exposure time when selecting USER MODE. **[0100]**

GAMMA: Setting gamma. **[MANUAL]**

Operation

PAGE 2 Setting tone of the image

- WB MODE Setting white balance. **[AWB]**
- AWB: White balance is controlled automatically.
- FULL OPEN: White balance is controlled without depending upon the condition of the subject.
- AWB HOLD: The existing white balance is held.
- USER MODE: Setting white balance manually. Adjust the value of USER R/G and USER B/G.
- USER R/G: Adjust the value of R/G when selecting USER MODE. **[02453]**
- USER B/G: Adjust the value of B/G when selecting USER MODE. **[07295]**
- OFFSET: Setting offset of the picture. **[000]**
- CONTRAST: Setting contrast of the picture. **[128]**
- SHARPNESS: Setting sharpness of the picture. **[H00 V00 COR00]**

PAGE 3 Setting the line marker

- GRAPHICS: Select whether the line markers are on or off. **[ON]**
- LINE: Select whether the line markers are displayed or not. **[ON]**
- LINE1 H: Adjust the horizontal line of the line marker 1; position by POS, width by SIZE and color by COLOR.
- V: Adjust the vertical line of the line marker 1; similarly.
- LINE2 H: Adjust the horizontal line of the line marker 2.
- V: Adjust the vertical line of the line marker 2.

Lines are displayed when both "GRAPHICS" and "LINE" are on.
Lines are not displayed when "SIZE" is set to 0.
ON/OFF of "GRAPHICS" is linked with ON/OFF of "GRAPHICS" of SHADOW MASK on PAGE 4.

PAGE 4 Setting the shadow masks

- GRAPHICS: Select whether the shadow masks are on or off. **[ON]**
- SHADOW: Select whether the shadow masks are displayed or not. **[ON]**
- GRADE: Adjust the grade of shading.
- H: Adjust horizontal masking quantity.
T: from the top, B: from the bottom
- V: Adjust vertical masking quantity.
L: from the left, R: from the right

Shadow masks are displayed when both "GRAPHICS" and "LINE" are on.
ON/OFF of "GRAPHICS" is linked with ON/OFF of "GRAPHICS" of LINE MARKER on PAGE 4.

Operation

PAGE 5 Setting other items

- RES/FPS: Setting resolution and frequency of output images.
Adjust corresponding your monitor screen. **[1080P 60]**

Select from formats below;

1080P 60, 1080P 30, 1080P 50, 1080P 25, 720P 60,
720P 50, 1080P 59.94, 1080P 29.97, 720P 59.94

- OSD SIZE: Changing size of the letters on the menu. **[LARGE]**
- PROFILE: Download the saved settings from the camera. **[PRESET 0]**
See "How to save" on the bottom of this page for saving.
- PATTERNS: Outputting test patterns. **[OFF]**
- GRAY: A gray scale chart is output.
- COLOR: A color pattern is output.
- GRAY+COLOR: A color pattern in the upper half of the screen and a gray scale chart in the lower half.
- IMAGE OUTPUT: Inverting images. **[STANDARD]**
- STANDARD: Not inverted images are output.
- H INVERSION: Horizontally inverted images are output.
- V INVERSION: Vertically inverted images are output.
- HV INVERSION: Horizontally and vertically inverted images are output.

PAGE 6 saving the settings

- EEPROM: Save the settings of PAGE1 to 5 in the camera.

How to save

Prior to saving, select the saving place (PRESET 0 to 7) of PROFILE on PAGE 5 and end setting on that PRESET *.
After various settings, move the cursor to SAVE of EEPROM on PAGE 6.
Depress the button D to display "Are You OK?" Depress the button D again to save. If depress the button C or E during a confirmation message is displayed, the settings are not saved.

* All the PRESETs are set to the initial values by the manufacturer.
Saving means overwriting them. Keeping one PRESET among PRESET 0 to 7 without any modification is recommended.

HOZAN TOOL INDUSTRIAL CO.,LTD.

1-2-12 Saiwaicho, Naniwa-ku, Osaka 556-0021, Japan
Tel : +81-6-6567-3111 Fax : +81-6-6562-0024