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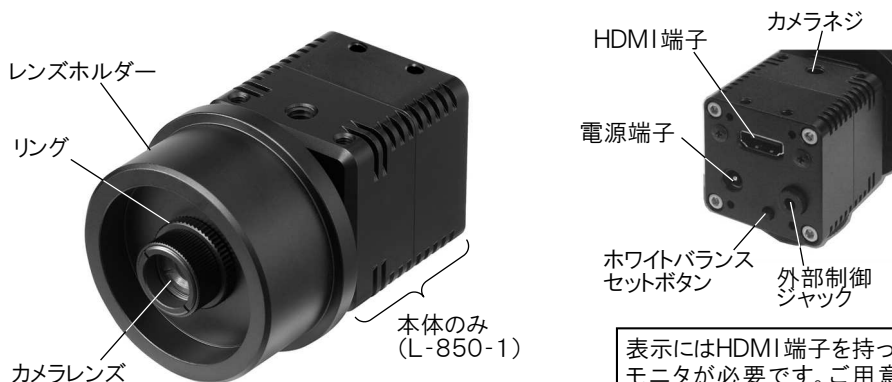
L-850 FULL HD CAMERA

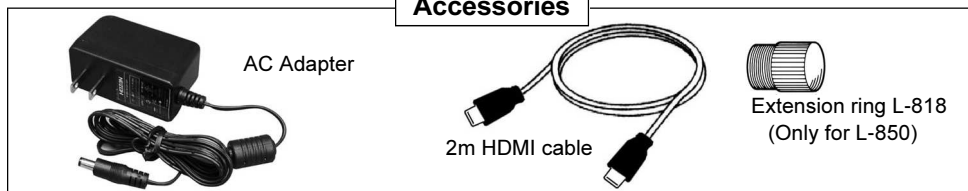
L-850-1 FULL HD CAMERA WITHOUT LENS

Thank you for purchasing the HOZAN L-850/L-850-1 FULL HD CAMERA. With proper care and handling, this fine instrument will provide years of trouble-free operation. Please read this entire instruction manual carefully before attempting to place this instrument in service. Please keep this instruction manual available for reference.

Identification of parts and contents

Accessories

Specifications

Element	1/2.8" 2.3 megapixel CMOS	Rating	12 Vdc
Image output effective pixels	1920(H) × 1080(V)	Power consumption	4.4 W
Video output	Based on DVI 1.0 RGB, 1080P60	Weight	228 g (without accessories)
Output connector	HDMI connector		
Lens mount type	C mount		

Warning and caution symbols

These symbols are used throughout the instruction manual to alert the user to potential safety hazards as follows :

⚠ Warning ... Notice when incorrect handling could cause the user's death or serious injury.

⚠ Caution ... Notice when incorrect handling could cause injury to the user or material damage.

Even if the instructions do not have **⚠ Caution** mark, there are some possibilities for a serious situation. Follow the instructions.

Precautions

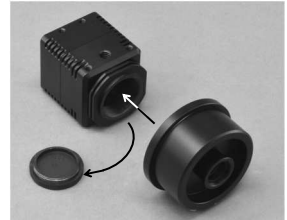
⚠ Caution

1. Use only the power adapter supplied with the camera, not any other alternative 12V source.
2. Do not modify the equipment in any way.
3. Do not expose the camera to moisture.
4. Avoid physical shock to the camera. This could cause failure to camera.
5. Do not expose the equipment to excessive heat, for example direct sunlight or other heat source.
6. Do not expose the lens of this camera to direct sunlight or other powerful light. Such exposure could damage the CMOS.

Preparation

Remove the cap of the camera, then screw the lens into the mount of the camera clockwise.

The connecting format of the camera and the lens is C mount. Using the L-815 ZOOM LENS with this camera makes precisely enlarged images. Also other C mount lenses in the market can be used.

Operation

- 1 Set up the L-850 on the appropriate equipment e.g. the L-521 STANDARD TABLE accompanying the L-509 HOLDER.

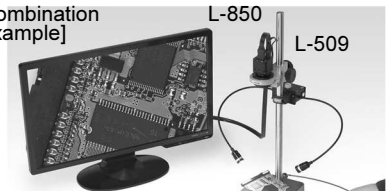
Also be connected to tripods and such in the market which have camera screws.

Camera screw

- 2 Install the L-850 on the holder or other equipment.

[Combination example]

L-521 STANDARD TABLE
The post is L-520 LONG POST.

Operation

- 3** Connect the unit and the monitor display, aspect ratio 16:9, which have HDMI inlet using the HDMI cable.

If use other than the provided HDMI cable

Double shielded HDMI cables specified as HDMI 1.3 or more and up to 5m length are recommended. The signal carried through a long cable beyond 5m may decrease and the monitor cannot display the image, depending also on the monitor's specifications. Please make sure of the standards and specifications of the cable and apparatus.

- 4** Insert the plug of the ac adapter into the power connector.
Select the input mode of the monitor side "HDMI".

[Regarding the magnification]

There is a ring, effective length 3mm, at the foundation of the optical lens. And the L-818, effective length 8.5mm, is accompanied (L-850 only). Various magnifications are gotten removing/attaching these rings. The relation between the combination and magnification is follows.

Combination	Magnification(×) *	Field of view mm	Working distance mm	Resolution at center** μ m
Optical lens singly (without ring)	9.4	54.5×30.5	330	88.5
Optical lens with ring	19	27.0×15.2	175	44.2
Optical lens with ring + L-818	46.5	11.1×6.2	85	17.5

* Using 24-type widescreen monitor. There is some variation depending on the characteristics of the TV to be used.

** The limit value at the center of field possible to recognize lines as lines which are drawn in the same thickness at the same interval.

- 5** While watching the monitor, focus above the object for best picture. When using the L-509, use the focus knob for precise focus after temporary focus by holder's up-down.

- 6** Remove the ac adapter from the outlet after use.

[Regarding white balance]

The white balance functions are altered among "full automatic" and "semi automatic" depressing the white balance set button longer than 2 seconds.

Full automatic function: The white balance is adjusted always automatically according to the photographing image.

Semi automatic function: The white balance is adjusted according to the image of the point in time when the white balance set button is depressed for an instant. This adjustment is held whether the image changes or not until the button is depressed again.

Using the optional L-815 camera controller, various settings can be done farther more. See page 4.

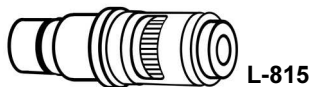
White balance
set button

Options

L-818 Extension ring
L-819 Extension ring

L-815 Zoom lens

If more large magnification is required, use extension rings singly or combining. The magnification, field of view and working distance are as follows.

Combination	Magnification(×) *	Field of view mm	Working distance mm	Resolution at center μ m
Optical lens (with ring) +L-818+L-818	75	7.0×3.9	61	11.0
Optical lens (with ring) +L-819	95	5.5×3.1	54	9.9
Optical lens (with ring) +L-818+L-819	122	4.3×2.4	45	7.8
Optical lens (with ring) +L-819+L-819	170	3.1×1.7	40	6.2
Optical lens (with ring) +L-818+L-819+L-819	195	2.6×1.5	38	5.5
L-815	27 to 185	19.4×10.8 to 2.8×1.6	105	31.2 to 4.9

* Using 24-inch widescreen monitor. There is some variation depending on the characteristics of the TV to be used.

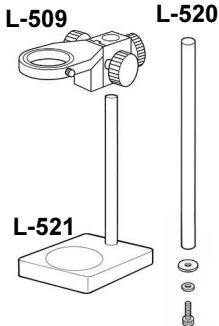
When use the L-818/819, setting a lighting system to compensate for reduced illumination caused by increased magnification is recommended.

L-509 Holder

L-521 Standard table

L-520 Long post

When inspect in low magnifying power(×9.4), exchange the post to the L-520 Long post.

L-850-2 Camera controller

Controls white balance, contrasts and others.

HOZAN TOOL INDUSTRIAL CO.,LTD.

1-2-12 Saiwaicho, Naniwa-ku, Osaka 556-0021, Japan
 Tel : +81-6-6567-3111 Fax : +81-6-6562-0024