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
L-498 EYEPIECE with SCALE

Thank you for purchasing the HOZAN L-498 EYEPIECE with SCALE. With proper care and handling, this fine instrument will provide years of trouble-free operation. Please read this entire instruction manual carefully before attempting to place this instrument in service. Please keep this instruction manual available for reference.


Identification of parts and contents

Dimensions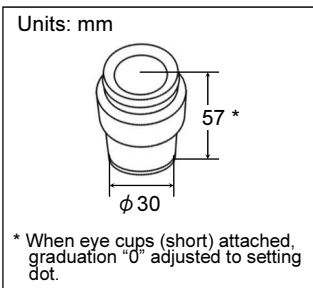

Micrometer, built-in

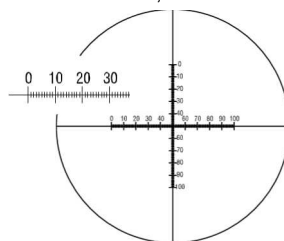(10 mm scale in 1/10 mm units)

Specifications

Magnification	10X
Field number	20
Weight	85 g with eyecup (short) attached
Applicability	L-46, L-51, L-461, L-514, L-KITs which include L-461 or L-514

Warning and caution symbols

These symbols are used throughout the instruction manual to alert the user to potential safety hazards as follows :

⚠ Warning ··· Notice when incorrect handling could cause the user's death or serious injury.

⚠ Caution ··· Notice when incorrect handling could cause injury to the user or material damage.

Even if the instructions do not have **⚠ Caution** mark, there are some possibilities for a serious situation. Follow the instructions.

Precautions


⚠ Caution

1. Do not disassemble or modify the eyepieces yourself, otherwise this could cause a malfunction.
2. Protect the unit against physical shock or vibration during transport. Otherwise, this could cause a malfunction.
3. Do not expose the unit to prolonged direct sunlight, high temperature, high humidity, or dust. Such exposure could cause a malfunction.

Preparation

- 1 First, make sure that the contents are complete.
- 2 Prepare an appropriate microscope (eyepieces of 10x are attached).
- 3 Exchange the existing eyepiece, left or right, with this eyepiece with scale, after loosen the eyepiece set screw.


Do not secure the eyepiece set screw yet.
- 4 Turn the eyesight revision ring to get a clear image of the micrometer.
- 5 Adjust focus according to the instruction manual of each microscope.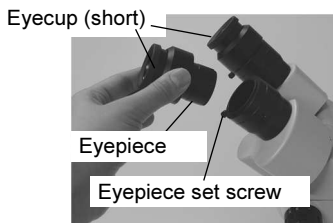

* Please refer to each instruction manual for operation of the microscope. Instruction manuals are available on our webpage. Utilize if lost.

Operation

- 6** Turn the eyepiece (micrometer) by its barrel according to the object if necessary.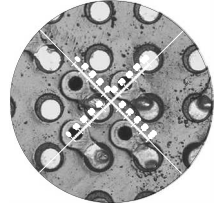

- 7** Secure the eyepiece set screw.

• Hints

The scale of this eyepiece is designed that a division is 0.1 mm when the magnification of the objective is 1x.

When the total magnification is 20x, on L-51 Stereo microscope/L-514 Stereo head scope, please twice the value.


On L-46 Stereo microscope/L-461 Stereo head scope, it can be used in the whole zoom area converting the value.

Please correspond a division to the real dimension looking a basic object such as a ruler if use for strict measurement.

Please use eyecups (long) if necessary.

< How to attach the eyecup (long) >

1. Remove the eyecup (short) .
2. Attach the eyecup (long) as shown in the illustration.

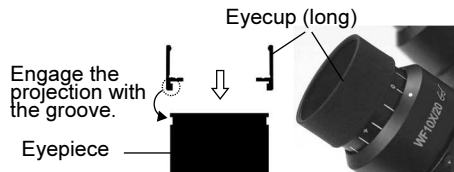Choose to suit

Eyecup (short)

- for persons in spectacles

Eyecup (long)

- for persons in contact lenses or without spectacles

Bringing inspector's cheekbones at eyecups (long) stabilizes his/her eye points.

- for blocking the slanting rays

Daily care

Lenses

- Wipe away any dust gently with a soft cloth or a piece of gauze.
- Clean away fingerprints or oil stains with a cloth which is lightly soaked in IPA or denatured alcohol (such as ethyl alcohol or methyl alcohol).
 - * Alcohol is extremely flammable, so handle it with care.

Note

- Do not allow any dust or fingerprints to get on the lens to get clear images.

Related products

Z-292

Antistatic cleaner

Static elimination of office automation and electrical appliances is tainted off.
Plastics and painted surfaces aren't infringed.
Prevents dust sticking to the lens by electrostatic influence.

Z-268

Rubber blower with compressor

The compressor strong/gentle can be selected.
Features an air inlet valve that prevents blowback from the inlet.
An accessory extension nozzle permits access to hard-to-reach places. Most suitable for microscopes which have lenses in cylindrical parts.

HOZAN TOOL INDUSTRIAL CO.,LTD.

1-2-12 Saiwaicho, Naniwa-ku, Osaka 556-0021, Japan
Tel : +81-6-6567-3111 Fax : +81-6-6562-0024