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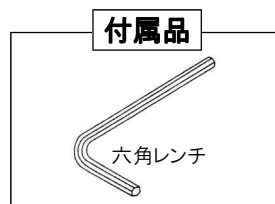
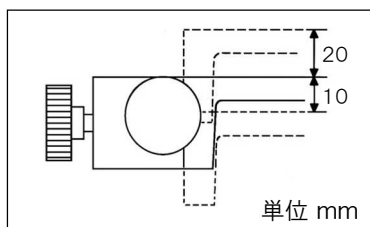
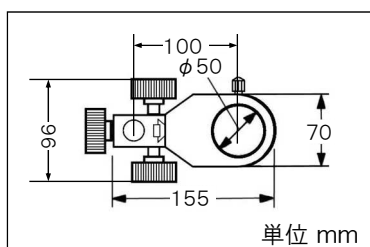
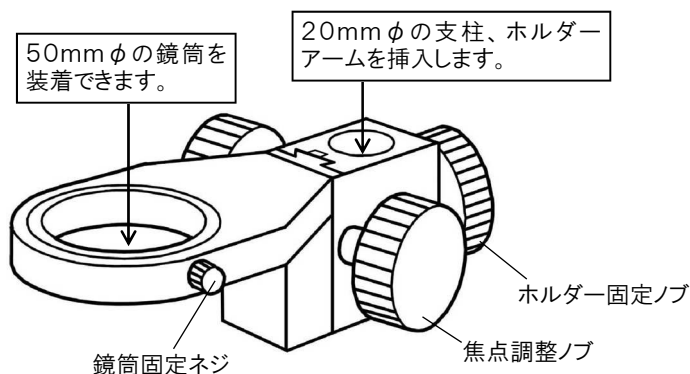
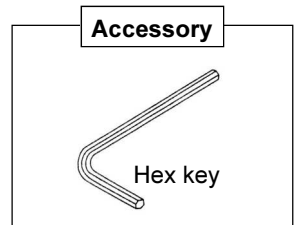
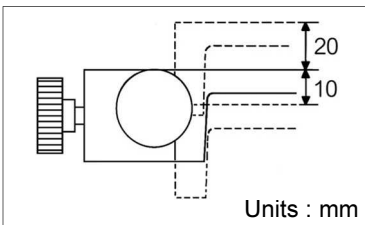
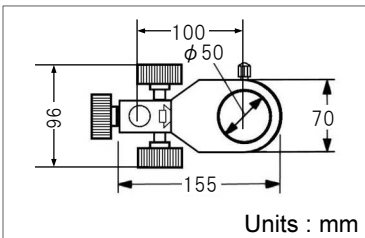
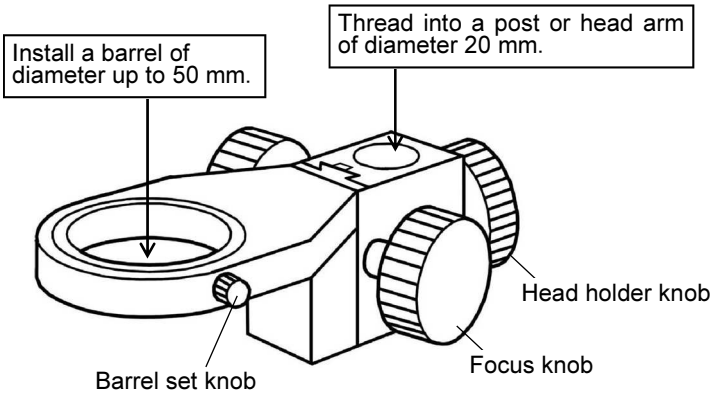
TEL(06)6567-3111 FAX(06)6562-0024

L-509 STEREO HEAD HOLDER

Thank you for purchasing the HOZAN L-509 STEREO HEAD HOLDER. With proper care and handling, this instrument will provide years of trouble-free operation. Please read this entire instruction manual carefully before attempting to place this instrument in service. Please keep this instruction manual available for reference.

Identification of parts and contents

Specifications

Dimensions	70(W)X50(H)X177(H)mm
Vertical stroke	20mm above 10mm below

Warning and caution symbols

These symbols are used throughout the instruction manual to alert the user to potential safety hazards as follows :

⚠Warning ... Notice when incorrect handling could cause the user's death or serious injury.

⚠Caution ... Notice when incorrect handling could cause injury to the user or material damage.

Even if the instructions do not have **⚠Caution** mark, there are some possibilities for a serious situation. Follow the instructions.

Precautions

⚠ Caution

Be sure that the head holder knob and barrel set knob are securely tightened before use, otherwise the head holder could slide down and be broken.

Operation

Prepare a base which is accompanied with a post of 20 mm dia.

Insert the L-509 into the post, and then secure by tightening the head holder knob at the position where the appropriate working distance can be gotten.

Precise focusing can be done with right and left focus knobs. That vertical stroke is 20 mm above and 10 mm below from the basic setting, the holder's top face and the block's top face are in the identical plane.

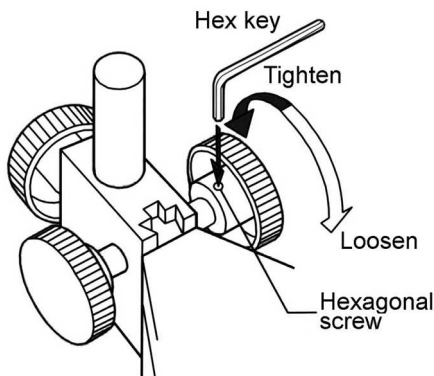
- * Refer to each instruction manual for the appropriate working distance, detailed uses of the microscope (head and eyepieces), camera and so on.

Adjusting tightness of the focus knob

The tightness of the focus knob for vertical movement can be adjusted.

Use the hexagonal key to loosen the screw at the neck of the focus knob, and then turn the knob clockwise to increase the tightness, or turn it counterclockwise to decrease the tightness. After adjusting, tighten the hexagonal screw.

- * Do not loosen unnecessarily, otherwise the stereo head will gradually slide down, which will interfere with your work.

HOZAN TOOL INDUSTRIAL CO.,LTD.

1-2-12 Saiwaicho, Naniwa-ku, Osaka 556-0021, Japan
Tel : +81-6-6567-3111 Fax : +81-6-6562-0024