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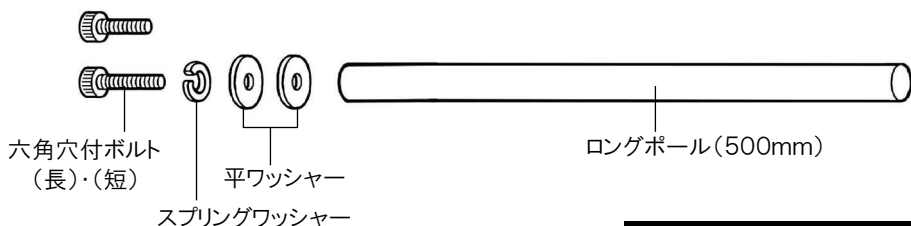
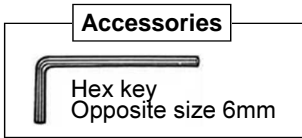
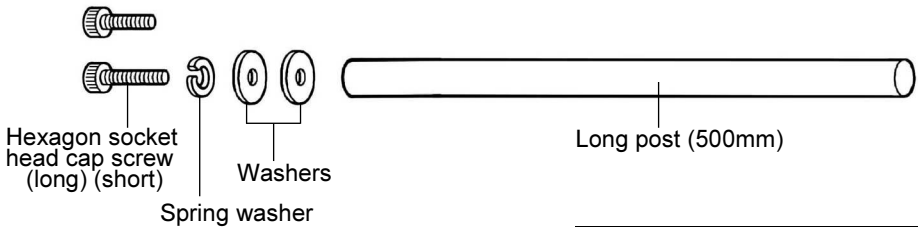
TEL(06)6567-3111 FAX(06)6562-0024

# L-520 LONG POST

Thank you for purchasing the HOZAN L-520 LONG POST. With proper care and handling, this instrument will provide years of trouble-free operation. Please read this entire instruction manual carefully before attempting to place this instrument in service. Please keep this instruction manual available for reference.

## Identification of parts

## Specifications

Dimensions	20mm $\phi$ $\times$ 500mm
Weight	1.3kg

## Warning and caution symbols

These symbols are used throughout the instruction manual to alert the user to potential safety hazards as follows :

**Warning** ... Notice when incorrect handling could cause the user's death or serious injury.

**Caution** ... Notice when incorrect handling could cause injury to the user or material damage.

Even if the instructions do not have **Caution** mark, there are some possibilities for a serious situation. Follow the instructions.

## Precautions

This post is for L-521 Standard table and L-525 Large table. Do not use for any other purpose.

**Caution**

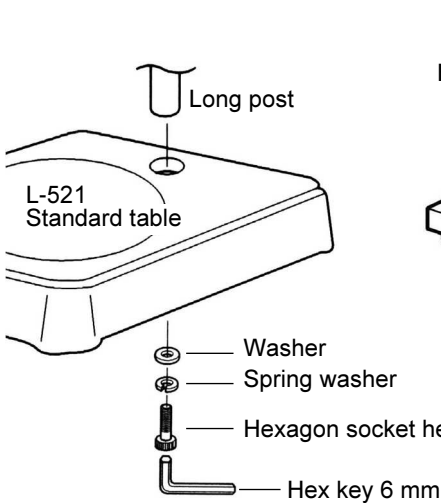
Be sure all parts are tightened securely, otherwise it could fall and the camera, microscope and the specimen could be damaged.

## Preparation for use


Prepare the L-521 STANDARD TABLE or L-525 LARGE TABLE for fixing this pole.  
Use the provided 6mm hex key.

- 1 Thread the hexagon socket head cap screw with the washer into the post fixing hole of the base upward from the bottom.  
Then join the post and secure the screw tightly.

[For L-521 Standard table]

[For L-525 Large table]

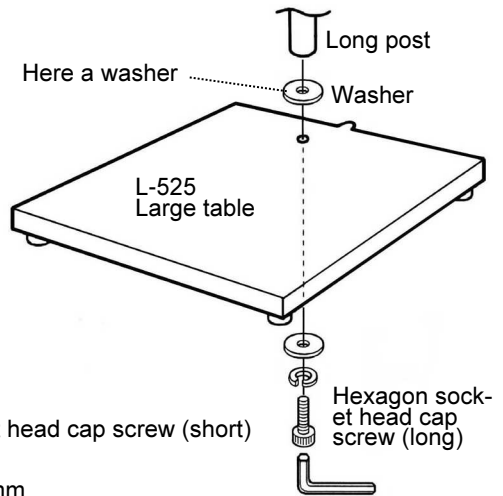- 2 Attach holders such as L-509 and L-804 on the pole to use cameras and microscopes.

### Caution

This is easily toppled by its height. Handle with care.

**HOZAN TOOL INDUSTRIAL CO.,LTD.**

1-2-12 Saiwaicho, Naniwa-ku, Osaka 556-0021, Japan  
Tel: +81-6-6567-3111 Fax: +81-6-6562-0024