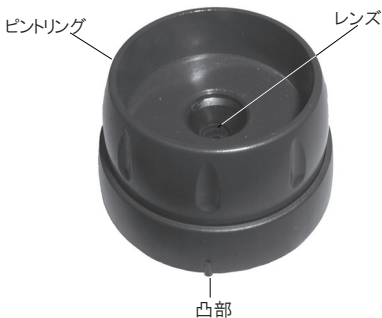
L-820-2

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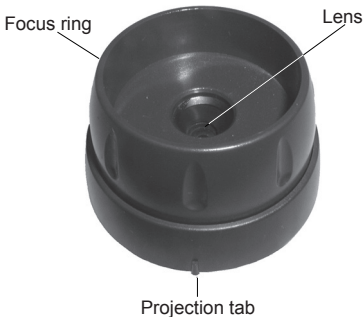
L-820-2

LENS UNIT

Thank you for purchasing the HOZAN L-820-2 LENS UNIT. With proper care and handling this fine instrument will provide years of trouble-free operation. Please read this entire instruction manual carefully before attempting to place this instrument in service. Please keep this instruction manual available for reference.


Identification of parts

Warning and caution symbols

These symbols are used throughout the instruction manual to alert the user to potential safety hazards as follows :

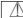 **Warning** ... Notice when incorrect handling could cause the user's death or serious injury.

 **Caution** ... Notice when incorrect handling could cause injury to the user or material damage.

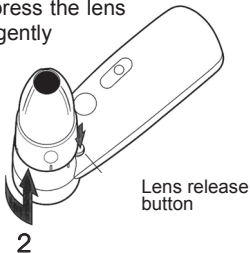
Even if the instructions do not have  **Caution** mark, there are some possibilities for a serious situation. Follow the instructions.

Precautions

This lens unit is for replacing the standard lens unit of the L-820 to acquire a normal (i.e., non magnified) picture. Do not use for any other purpose.


Operation

- 1 To remove the standard lens unit from the L-820, depress the lens release button and gently turn the lens counterclockwise while pulling.


Operation

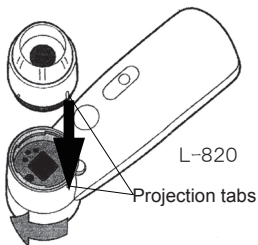- 2** To connect this lens unit to the L-820, bring the projection tabs on both the lens unit and the L-820 into alignment. Gently insert the lens unit into the L-820, then turn it clockwise until it locks into position.

- 3** You can adjust the focus by rotating the focus ring.

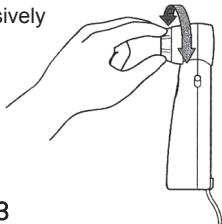Be sure not to excessively turn the focus ring in a counterclockwise direction, as it may fall out of the unit.

Operation

The illumination setting for the application software can be used just as it is set for the standard lens unit. However, make any fine adjustments you think are necessary.

For operating the application software of the L-820, choose the "On Screen Manual" tab from the menu in the CD-ROM that was provided with the L-820.

HOZAN TOOL INDUSTRIAL CO.,LTD.

HEAD OFFICE : 1-2-12 Saiwaicho, Naniwa-ku, Osaka 556-0021, Japan

Tel : (06)6562-0023

Fax : (06)6567-3115