L-845

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L-845 ADAPTER KIT

Thank you for purchasing the HOZAN L-845 ADAPTER KIT. With proper care and handling, this fine instrument will provide years of trouble-free operation. Please read this entire instruction manual carefully before attempting to place this instrument in service. Please keep this instruction manual available for reference.

Identification of parts and contents

Dia. of eyepiece	Combination
φ 23.2 mm	①
φ 30 mm	①+②
φ 30.5 mm	①+③

This is a lens system for adapting cameras which have C mounts to HOZAN's microscopes. Combine corresponding diameter of the eyepiece as the above table.

Warning and caution symbols

These symbols are used throughout the instruction manual to alert the user to potential safety hazards as follows :

⚠ Warning ... Notice when incorrect handling could cause the user's death or serious injury.

⚠ Caution ... Notice when incorrect handling could cause injury to the user or material damage.

Even if the instructions do not have **⚠ Caution** mark, there are some possibilities for a serious situation. Follow the instructions.

Precautions


⚠ Caution

1. Do not disassemble or modify the products. This could cause malfunction.
2. Avoid water.
3. Handle with care. Do not drop down. Avoid physical shock.

How to attach to the microscope

- 1 Recognize the eyepiece's outer diameter of the microscope to be used. Remove one of eyepieces and actually measure its part which is snugged into the microscope.

Since this device will be inserted to the barrel of the microscope, that removed eyepiece is not used. Store separately.

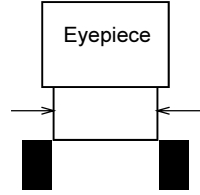- 2 When the measurement result is 30 mm or 30.5 mm, attach the corresponding conversion tube each.

If the measurement result is 23.2 mm, skip over this procedure.

- 3 Attach to the USB camera, screwing clockwise until snug.

- 4 Insert into the left barrel of the microscope.

Caution

Using this device makes the focal length different. Therefore refocusing on every magnification is necessary.

Combination example
L-802+L-845+L-51
+L-711+L-527

HOZAN TOOL INDUSTRIAL CO.,LTD.

1-2-12 Saiwaicho, Naniwa-ku, Osaka 556-0021, Japan
Tel : 81-6-6567-3111 Fax : 81-6-6562-0024